



Customer Guidance: Windows Settings to Maximize Communication Efficiency with SEAL Instrumentation

Instructions:

- Open the Control Panel
- Click on Power Options. If Power Options isn't an immediately available option, navigate to System and Security, then Power Options.
- Perform the following steps for both Balanced and Power Saver
 - Click Change plan settings
 - Click Change advanced power settings
 - Expand the Hard disk option
 - Expand the Turn off hard disk after and set to Never (0 minutes)
 - Expand the Sleep option
 - Set Sleep after and Hibernate after to Never (0 minutes)
 - Expand the USB Settings
 - Set USB selective suspend setting to Disabled
 - Click OK in the Advanced Power Settings window
 - Click Save changes in the Edit Plan Settings window
 - (Perform these steps for both Balanced and Power Saver)
- Navigate back to the Control Panel menu
- Click on Folder Options. If Folder Options isn't an immediately available option, navigate to Appearance and Personalization, then Folder Options
- Select the View tab
- Deselect the Use Sharing Wizard (Recommended) option
- Click OK